


1. Take sample from a faucet that is not leaking. Remove any attachments from the faucet.



2. Run water at full force for 5 min. to flush out pipes. Turn off water and flame tap with a lighter or match to sterilize it.



3. Turn water back on, but not full force. Tear Whirlpak bag at perforation being careful to avoid contaminating the inside edge of the bag. Grasp white tab at side of bag and carefully open bag. Fill bag 4/5 full; then, holding YELLOW wire tabs, **whirl bag end over end 3 complete revolutions.**

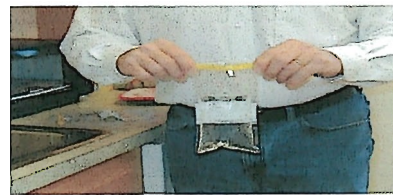

4. Turn tabs up **and twist them securely around each other.** Label bag.



5. **CHILL SAMPLE**, make sure lab receives sample within 30 hours.